

Surgery tables, Heating and Gas evacuation systems for mice and rats for Veterinary use

Surgery tables

Controlled heating systems

Gas evacuation systems

Manageable

Easy to use

combi-vet®



Rothacher Medical GmbH



SURGERY TABLES, HEATING- AND GAS EVACUATION SYSTEMS FOR

Overview


Especially designed for research field, we have different surgery tables, Heating- and anesthesia gas evacuation systems available. With these different accessories you will find a perfect solution to perform your surgeries. These systems can be added to existing or to our combi-vet anesthesia system.

combi-vet® Surgery table standard

For research, we have a standard surgery table with side holder for elastic rubbers or with a left or right side metal frame for FST clamps and holders. Please consult the FST General Catalogue. The table comes with a mask system for mice or small rats. Size 190 x 270 mm, working space 190 x 140 mm, 20 mm thick, easy washable as well as autoclavable. An optional fan with power supply can be connected to the downdraft anesthetic gas.



Order-Nr. PS-0451	Surgery table standard
Order-Nr. PS-0451-FST	Surgery table FST
Order-Nr. PS-0452	Fan with power supply 110-240 V/1,5 A DC

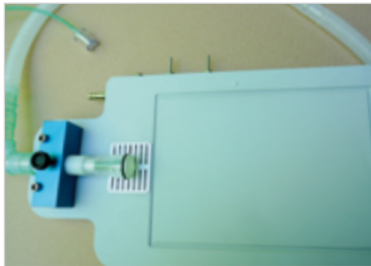

Order-Nr. PS-0451



Order-Nr. PS-0451-FST




Order-Nr. PS-0452


combi-vet® Surgery table standard with Physitemp temperature controller unit

The heating plate is controlled by a Physitemp temperature controller unit and a rectal probe is controlling the body temperature on mice or rats. The size of the heating plate fits our standard surgery table PS-0451 or PS-0451-FST

Order-Nr. TCAT-2LV	Temperature controller T-CAT
Order-Nr. HP-1M	Heating plate for T-CAT System
Order-Nr. RET-3 ISO	Isolated Rectal Probe for mice

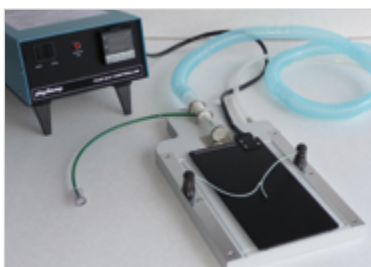

Order-Nr. TCAT-2LV

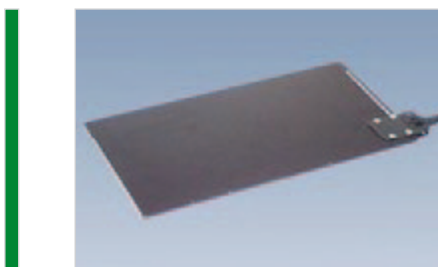

Order-Nr. HP-1M




Order-Nr. RET-3 ISO

R MICE AND RATS FOR VETERINARY USE

combi-vet® Surgery table with metallic frame and Safecontrol
«Economy» Temperature controller for mice

3 features - all in one Surgery table, Heating system and Gas evacuation

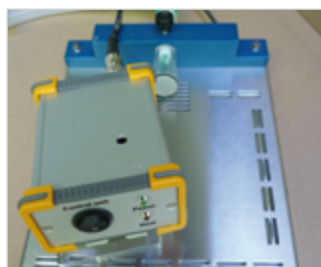

Another surgery table is our new developed set. The modular system can be step by step adjusted to your needs. It is available with a Temperature Controller and a incorporated heating pad build in under the metallic frame. It will deliver a constant temperature between 36 C - 38 C. The size is 190 x 230 mm and the working space is 100 x 100 mm, 20 mm thick. The stainless steel plate is 1,8 mm thick. A mask system for mice & small rats is mounted on a metal holder to the table. In addition, the anesthetic gas will downdraft from each side and under the mask. The exposure for users to anesthetic gas will be reduced to a very low level.

Module 1 Surgery table

Order-Nr. PS-0703 SM

Surgery table small for mice without heating plate


Order-Nr. PS-0703 HP

Surgery table small for mice with heating plate

Order-Nr. PS-0704

Metal holder mounted on plate with Mask for mice or small rats

Mask diameter is 20 mm , It will fit PS-0703 SM and HP version

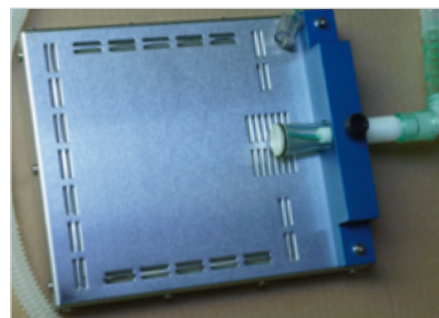

Module 2 Gas evacuation fan system

Order-Nr. PS-0452 Set

Fan with on/off switch on cable, without power supply



Module 3 Temperature controller

Order-Nr. PS-070-A

Temperature Controller unit with Power

Supply 110-220 V 1,7 A EN 60601/1



Order-Nr. PS-0545

Power supply only 110-240 V/1,7A EN60601/1



combi-vet® Surgery table with metallic frame and «Safecontrol» Temperature controller for mice & rats and reptiles

3 features - all in one Surgery table, Heating system and Gas evacuation

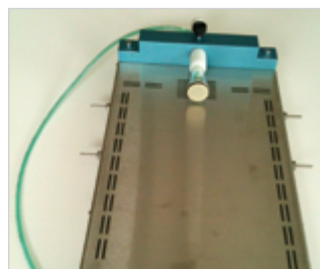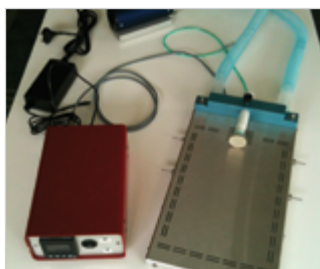

This surgery table is available as a modular system. Compared to the Safecontrol «economy» model for mice the system has a very accurate Temperature Controller, a larger surgery table and can be used for mice & rats or reptiles. The Safecontrol unit maintain and heat the surface between 35 C - 45 C very accurate. It can be adjusted very easily and a rectal probe is not necessary. The size of the surgery table is 240 x 190 mm with a working space of 170 x 110 mm. 18 mm thick. Enough space to fit bigger rodents. The stainless steel plate is 1,8 mm thick and easy to clean. A mask system for mice and rats is mounted on a metal holder to the surgery table. In addition, the anesthetic gas will downdraft from each side and under the mask. The exposure for users to anesthetic gas will be reduced to a very low level. The system makes the working area more safer for any user.

Module 1 Surgery table

Order-Nr. PS-0703 LG


Surgery table large without heating plate

Order-Nr. PS-0703 LG HP

Surgery table large with heating plate

Order-Nr. PS-0704 LG

Metal holder mounted on plate with Mask for mice or rats Mask, diameter is 20 mm or 30 mm exchangeable, It will fit PS-0703 LG and HP version

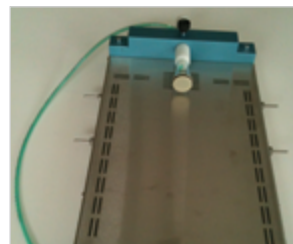

Module 2 Gas evacuation fan system

Order-Nr. PS-0452 Set

Fan without power supply with switch an/off on cable (Power supply source from Temperature Controller)



Module 3 Temperature controller

Order-Nr. PS-071-A


Temperature Controller unit programmable with Power Supply 110-220 V 24 V / DC 2,5 A with 12 V DC Outlet

Order-Nr. PS-0545

Power supply only 110-240 V/ 24 V DC 2,5 A



More accessories? Please request our 20 page documentation.

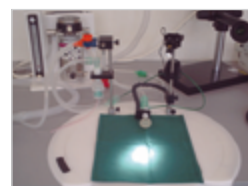

We reserve anytime the right modifications!